Referencia: PA0045

Ficha Técnica

Producto: **BRILLIANT GREEN AGAR (BGA)**

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Ficha Técnica

Producto: **BRILLIANT GREEN AGAR (BGA)**

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- ATLAS, R.M. & L.C.PARKS (1993) Handbook of Microbiological Media CRC Press. BocaRaton. Fla. USA.
- CLESCERI, L.S., A.E. GREENBERG & A.D. EATON (Eds) (1998) Standard Methods for the Examination of Water and Wastewater 20th ed. APHA- AWWA-WEF Washington DC. USA.
- DOWNES, F.P. & K. ITO (2001) Compendium of Methods for the Microbiological Examination of Foods.4th ed. APHA. Washington DC.
- EUROPEAN PHARMACOPOEIA (2006) 5th edition. Suppl. 5.6 EDQM. Council of Europe. Strasbourg. USA.
- FORBES, B.A., D.F SAHM & A.S. WEISSFELD (Eds) (1998) Bailey & Scott's Diagnostic Microbiology 10th ed. Mosby. St Louis, Mo. USA.
- HORWITZ, W. (2000). Official Methods of Analysis of the AOAC Internacional 17th ed. Gaithersburg. MD. USA.
- ISO/TS 11133-1: 2009. Microbiology of food and animal feeding stuffs.- Guidelines on preparation and production of culture media. Part 1: General guidelines on quality assurance for the preparation of culture media in the laboratory.
- ISO/TS 11133-2: 2003 Corr. 2004. Microbiology of food and animal feeding stuffs.- Guidelines on preparation and production of culture media. Part 2: Practical guidelines on performance testing of culture media.
- KAUFFMAN, F. (1935) Weitere Erfahrungen mit der kombinierten Anreicherung verfahren für *Salmonella* bazillen. Z. Hyg. Infekt. Krhn, 117; 26-32.
- MacFADDIN, J.F. (1985) Media for isolation-cultivation-identification- maintenance of medical bacteria. Williams & Wilkins. Baltimore. Md. USA.
- MURRAY, P.R., E.J. BARON, J.H. JORGENSEN, M.A. PFALLER & R.H. YOLKEN (Eds) (2003) Manual of Clinical Microbiology 8th ed. ASM Press. Washington DC, USA.
- OSBORN, W.W. and STOKES, J.L. (1955) The determination of Salmonellae in Foods. Ottawa: Food and Drug Laboratoires. 1962.
- US FDA (Food and Drug Administrations) (1998). Bacteriological Analytical Manual. 8th ed. AOAC International. Gaithersburg, MD. USA.